

Formalin Detection Test Kit in Food



Number of Test / Kit: 10 Test

Limit Detection: 0.5ppm

Shelf-life: 2 Years at room temperature

Procedure:

1. For liquid sample, pour the sample into the bottle of Formalin test reagent 1 upto 1/3 of the bottle. In case of solid sample, rinse the sample with sufficient clean water and transfer the rinsed water into the Formalin test reagent 1, close the cup and shake to completely dissolve the solid reagent.
2. Transfer the content from the bottle of Formalin test reagent 1 into the bottle of Formalin test reagent 2, close the cap and shake.
3. Transfer the content from the bottle of Formalin test reagent 2 into the bottle of Formalin test reagent 2, close the cap and gently swirl to mix the content, notice the color

Result Evaluation:

If the liquid turn pink or red, indicates that sample contains formalin.

Caution:

- Wash with running water and soap upon contact
- Keep the test kit out of children reach.